## Quick Start Op-DigiVac Scientific Measurement & Control

## Automated High Vacuum Pump-down System with SNAP Vacuum Control

## How to Install the Nitrogen/Inert Gas Tank:

- 1.On the back of the cart where the system connects, take the tape off the shelf that is folded up.
- 2. Place the tank (customer supplied, 40 ft³ recommended) on the rubber pad that is attached to the shelf.
- 3. Use the strap of the bracket around the tank and tighten it to secure the tank in place.
- 4. Attach the Regulator to the top of the tank. Make sure that the Regulator is set at no more than 1 psig.
- 5. Connect the regulator via hose to the KF16 fitting on the cart.



